



Biochemicals and Ready-to-Use Buffers

New!

Pre-mixed Buffers available in **Sachets**

- ✓ **Compact** for ease of storage
- ✓ **Huge savings** on shipping costs
- ✓ **Simple to Use**, just add de-ionised water to obtain 10X buffer
- ✓ **Suited for** Protein/Nucleic Acid Analysis, Electrophoresis and Blotting



Product Information

Product Number	Description
BUF-3009-10X1L	10X Tris-Acetate-EDTA (TAE) Buffer, Premixed Powder (sachet), Ultra Pure Grade, 1L
BUF-3019-10X1L	10X Tris-Borate-EDTA (TBE) Buffer, Premixed Powder (sachet), Ultra Pure Grade, 1L
BUF-2049-10X1L	10X Phosphate Buffered Saline (PBS), Premixed Powder (sachet), Ultra Pure Grade, 1L

Instruction for Use



Open sachet and empty entire content



Dissolve powder completely in 1L of distilled de-ionised water to yield a 10X buffer solution



Dilute buffer to desired concentration before use (autoclave/ filter if desired)

For ordering and enquiries, please contact

1st BASE Pte Ltd
41 Science Park Road #01-22/23/28
The Gemini, Singapore Science Park II
Singapore 117610

Tel: +65 6775 7318
Fax: +65 6775 7211
Email: info-sg@base-asia.com

First BASE Laboratories Sdn Bhd
No 7-1, Jalan SP 2/7, Tmn Serdang Perdana
Seksyen 2, Seri Kembangan 43300
Selangor Darul Ehsan, Malaysia

Tel: +603 8943 3252
Fax: +603 8943 3243
Email: info-my@base-asia.com